



***Kem Medical Products Corp.***

***KemSure<sup>®</sup> Formaldehyde  
Neutralizing Powder***

**Catalog # 9144**

**Directions:**

1. Pour **KemSure<sup>™</sup>** Formaldehyde Neutralizing Powder into container with waste formaldehyde solution which you intend to neutralize (one container neutralizing powder per ¼ gallon waste solution).
2. Stir solution to mix.
3. When reaction is complete (30 minutes), the neutralized solution is then safe for disposal.

*Use the entire contents of this container in one application. This product is not designed to be used in smaller portions.*

*Note:* Always maintain adequate ventilation in accordance with applicable regulations. Disposal is subject to federal, state, and municipal sewer regulations.

**Precaution:**

This product is not to be ingested or brought into contact with the skin and eyes. Review manufacturer's Safety Data Sheet. Keep away from children.